

Equipment Start Up and User Training Check List G 7881

Please complete and submit back to Miele Technical Support Department
 Fax # 800-220-1348 or email ProService@Mieleusa.com

Water Quality Test:

Hardness		Grains Per Gallon
Chlorides		Milligrams Per Liter

Installation Check - In

Inspected

1. Check the water lines are not kinked or bent to restrict flow	
2. Check in-line water filters are properly placed in stainless steel adapters	
3. Check water supply for Final Rinse in properly set (either tap or DI water)	
4. Check electrical supply is provided as specified	
5. Check drainage is provided as specified	
6. Check machine is level or secure on base if applicable	
7. Check all inserts are adjusted correctly - Top spray arm flange	
8. Check liquid detergent containers are full and operating properly	

Operator Training Check List

Please explain and demonstrate the following:

Performed

1. Control panel indicator lights and display	
2. Water softener electronically set and regeneration initiation	
3. Powder detergent door dispenser operation	
4. Liquid dispenser volume setting control - primed system	
5. Neutralizer dispenser volume setting control - primed system	
6. Mieleclear door container filled	
7. Triple circulation filters, removal and cleaning	
8. Review reset switch and minor troubleshooting procedures	
9. Review preventative maintenance procedures	
10. Perform test run with end-user(s)	
11. Explain use of baskets & inserts, adjustment of couplings	
12. Program selector dial and program cycle definitions	
13. Drying selection - explain drying operation	
14. Reviewed instrument suitability guidelines	

Personnel Training List: List all present during the training

Date of Start-Up:	Serial Number:
Client Name:	

For any additional questions please contact
Miele Professional: (800) 991-9380
9 Independence Way
Princeton, NJ 08540